


## Assembly of passive samplers for BTEX

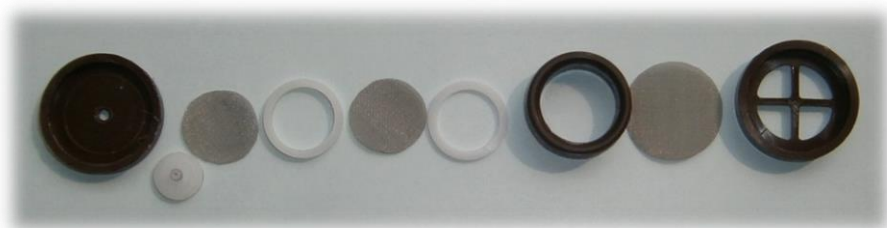

1. Wash all parts of passive samplers by tap waters. Rinse twice by deionized water. Leave on the air to dry. Do not dry by heat.

2. Attach the thick O-ring into the bottom of the body.



3. Place SS mesh on the O-ring.



4. Place thin O-ring on the SS mesh.



5. Join the body and the bottom.



6. Place the big SS mesh into the cap.



7. Join the cap and the body.



8. Fill PS with charcoal through the hole on the bottom.



9. Cap tightly. Store PS in aluminum zipper envelop.



- Wear glove when working with passive samplers and assembly should be done in clean air to avoid contamination.
- Store ready-in-use passive samplers in clean air at low temperature.
- Send passive samplers to Enviochem Binh lan for analyses, washing and replacement absorption medium.