



ENVIKIT pH

ENVIKIT 01pH  
4.0 – 9.5, 0.5 increment




- Thank you for choosing products of Binh Lan ENVIBIOCHEM. Please read instruction manual carefully before proceeding to ensure accurate results.
- Always cap reagent bottle gently to avoid evaporation.



1. Fill the tube with 3 ml (1/2 of the tube) of sample water.



2. Add 2 drops of the Indicator. Cap the tube and invert to mix.

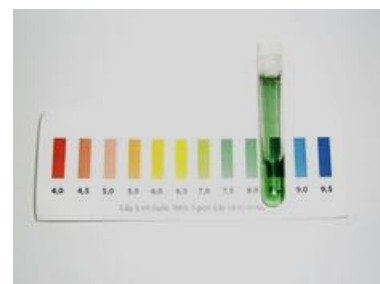

3. Find the best matched color in the color chart. Record pH.

## BINH LAN ENVIBIOCHEM COMPANY Limited

**Web:** <http://www.binhlan.com>      **Email:** [binhlanbiochem.gmail.com](mailto:binhlanbiochem.gmail.com)

**Bank account:** ACB Bank, Binh Lan Company, Account number: 1246888

**Office 1:** Science-Technology Business Incubator of HoChiMinh City  
268 Ly Thuong Kiet Street, District 10, HoChiMinh City, Vietnam

**Office 2:** 595/208A Cach mang thang 8, Ward 15, District 10, HoChiMinh, Vietnam

**Tel:** 08-39700106, **Hotline:** 0985714580

For technical support, quotation and order

Email to [binhlanbiochem.gmail.com](mailto:binhlanbiochem.gmail.com)